

# EPA Method TO-17 Sampling Instructions



1. Wear gloves whenever handling the sorbent tubes
2. Open container and remove foil from around the sorbent tube
  - a. Retain both for return shipping to the laboratory
3. Use a 9/16 inch and a 1/2 inch open ended wrench to remove caps
4. UN-Grooved end of the sorbent tube is attached to the pump – Arrow indicates direction of air/sample flow
5. Do not load with more than 1 to 4 Liters of sample
6. Once sampling is complete, remove sorbent tube from pump and recap sorbent tube
7. Wrap each sorbent tube in foil and place in original container
8. Ship on ice

If at any point you have questions, please call your Calscience Project Manager at 714.895.5494

*The difference is service*

